


# *Red*

## Strong Fluorescence Imager

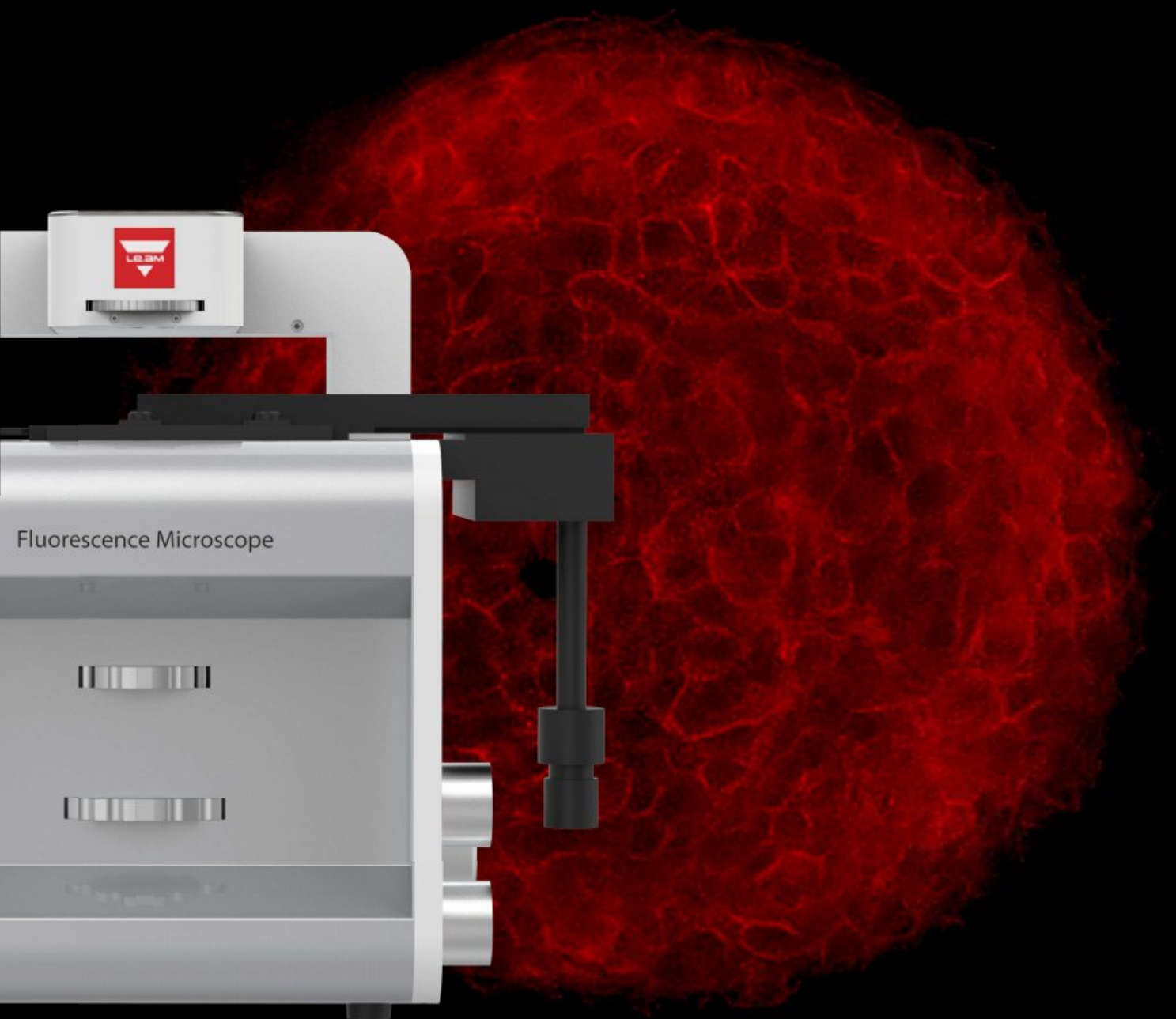



# Red

Strong Intensity  
Strong Signal  
Fluorescence




## Strong Optical Illumination Path


Strong fluorescent signal from optimized illumination path.  
Maximizes visualization of delicate biological samples.

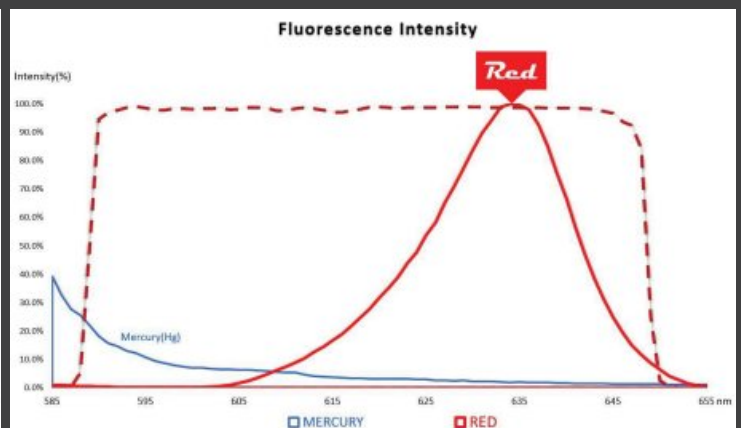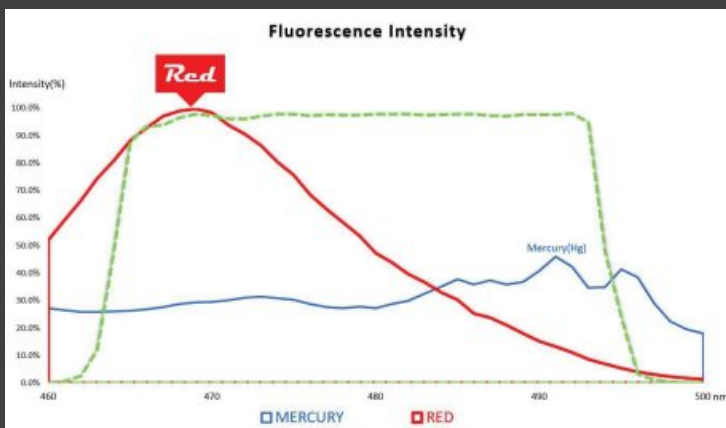

A strong optical illumination path results in strong intensity and produces a strong fluorescence signal

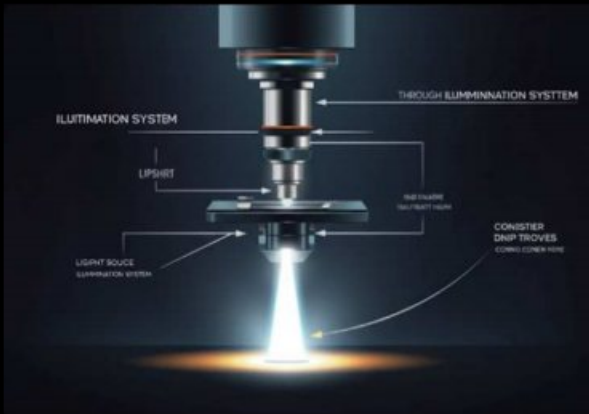

Fluorescence microscopy utilizes highly sensitive detectors to capture even the faintest of fluorescent signals, enabling the visualization of even the most subtle cellular features.

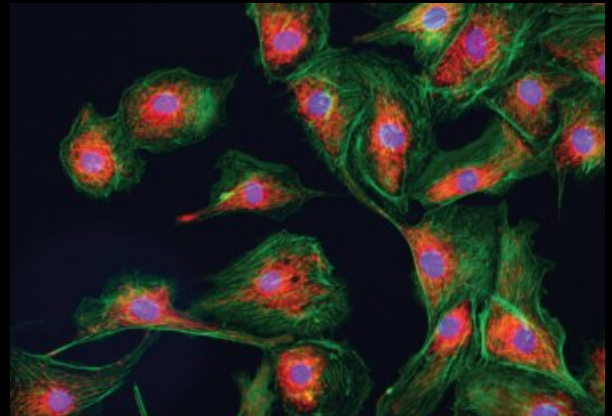


### Strong Signals


The use of fluorescent probes and markers provides a powerful amplification of target signals, allowing for the clear identification and tracking of specific biomolecules and structures.

Sometimes, you don't need a confocal microscope


Confocal

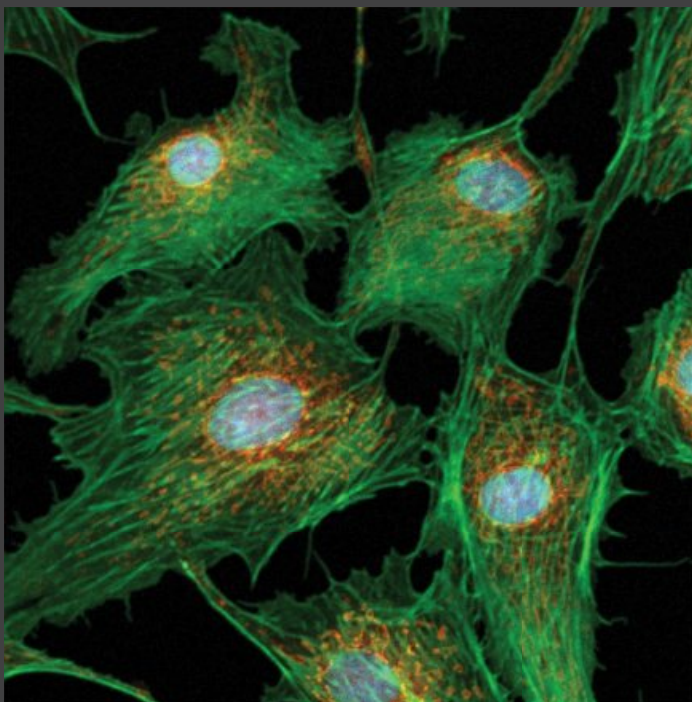

*Red*

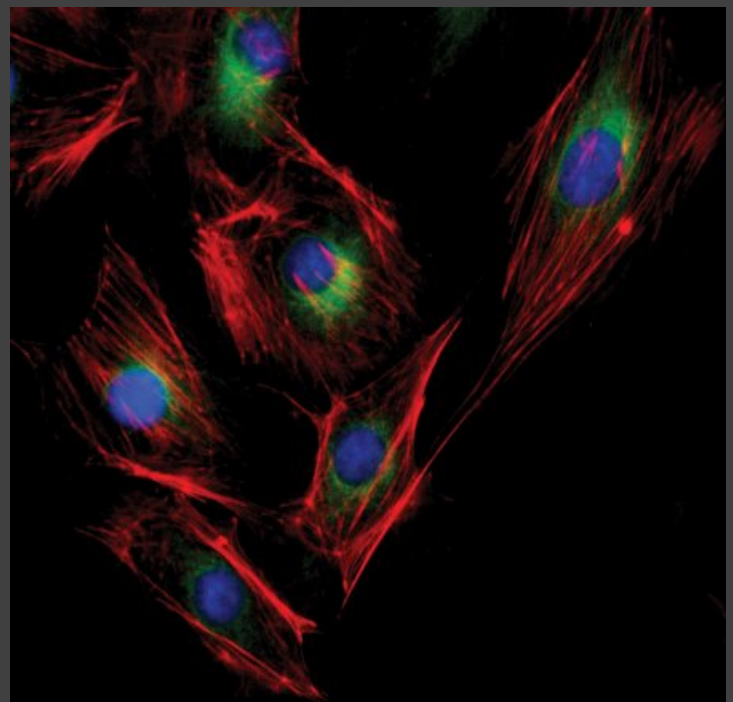



- Confocal microscope : FluoCells™ Prepared Slide #1 (Mito Tracker™ Red CMXRos / Alexa Fluor™ 488 Phalloidin/DAPI)
- *Red* Fluorescence imager : FluoCells™ Prepared Slide #2 (Mouse Anti- $\alpha$ -tubulin/AlexaFluor™ 488 Texas Red™-X Phalloidin/DAPI)

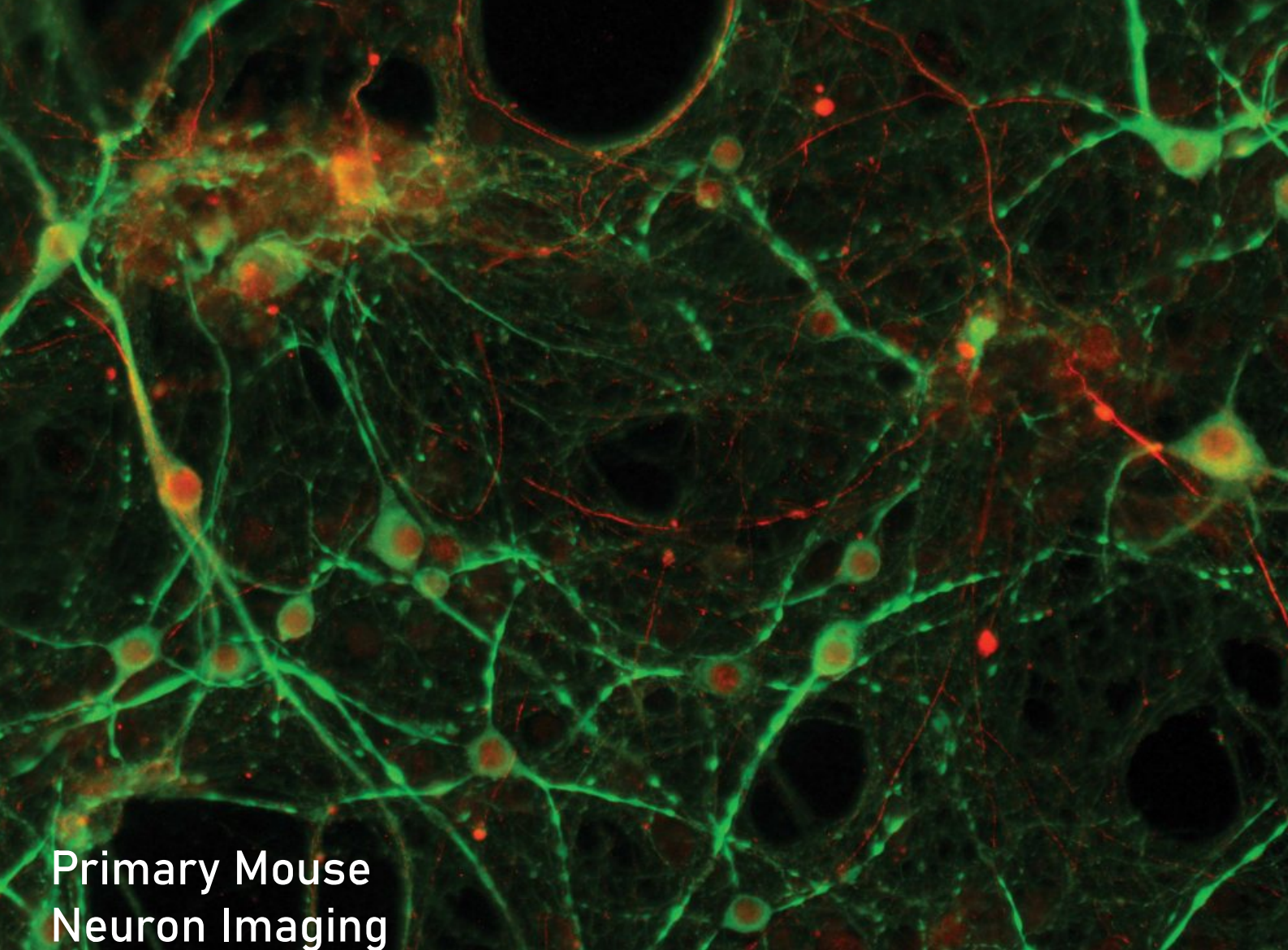

## Primary Mouse Neuron Imaging

Analyzing these fluorescent markers in primary neuron cultures helps elucidate the cellular mechanisms underlying neurodegenerative diseases like Parkinson's.

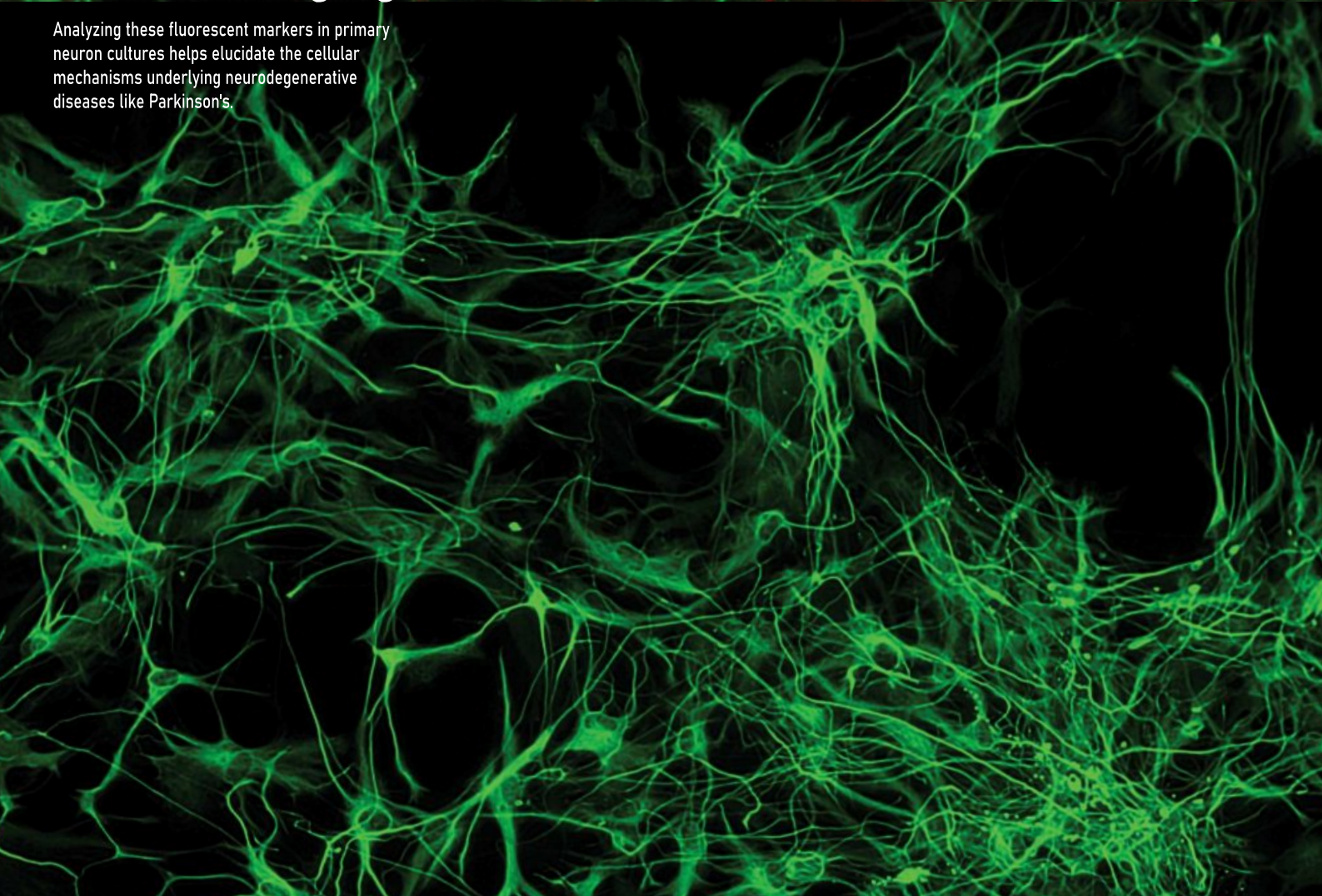

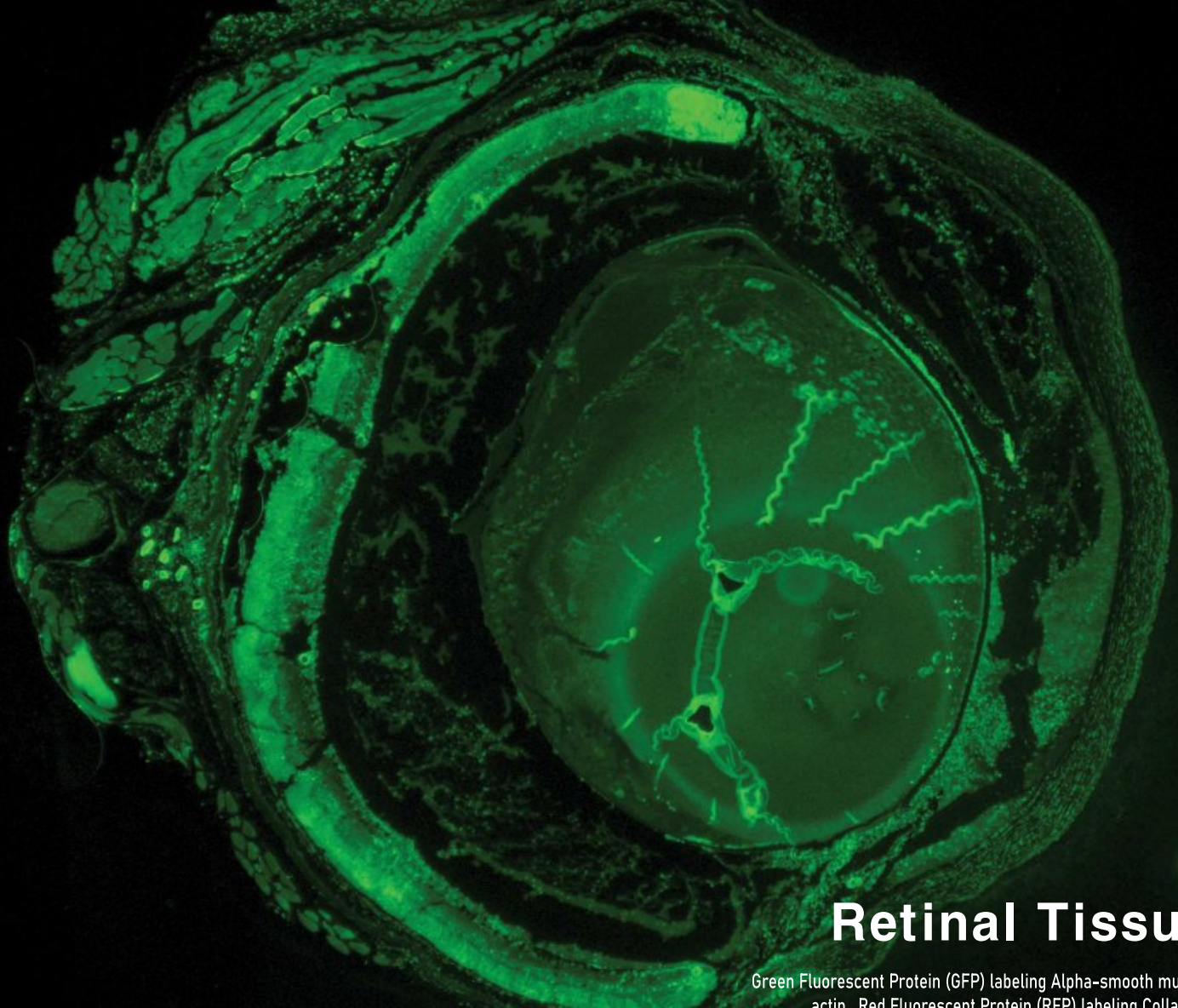



## Retinal Tissue

Green Fluorescent Protein (GFP) labeling Alpha-smooth muscle actin , Red Fluorescent Protein (RFP) labeling Collagen1  
4X objective for imaging Paraffin section of retina, 3 $\mu$ m thickness.

## Olfactory Bulb


The olfactory bulb is a crucial structure in the brain that processes and transmits smell information. This DAPI-stained image, captured using a 10X objective, image stitching, showcases the intricate cellular organization of this region.

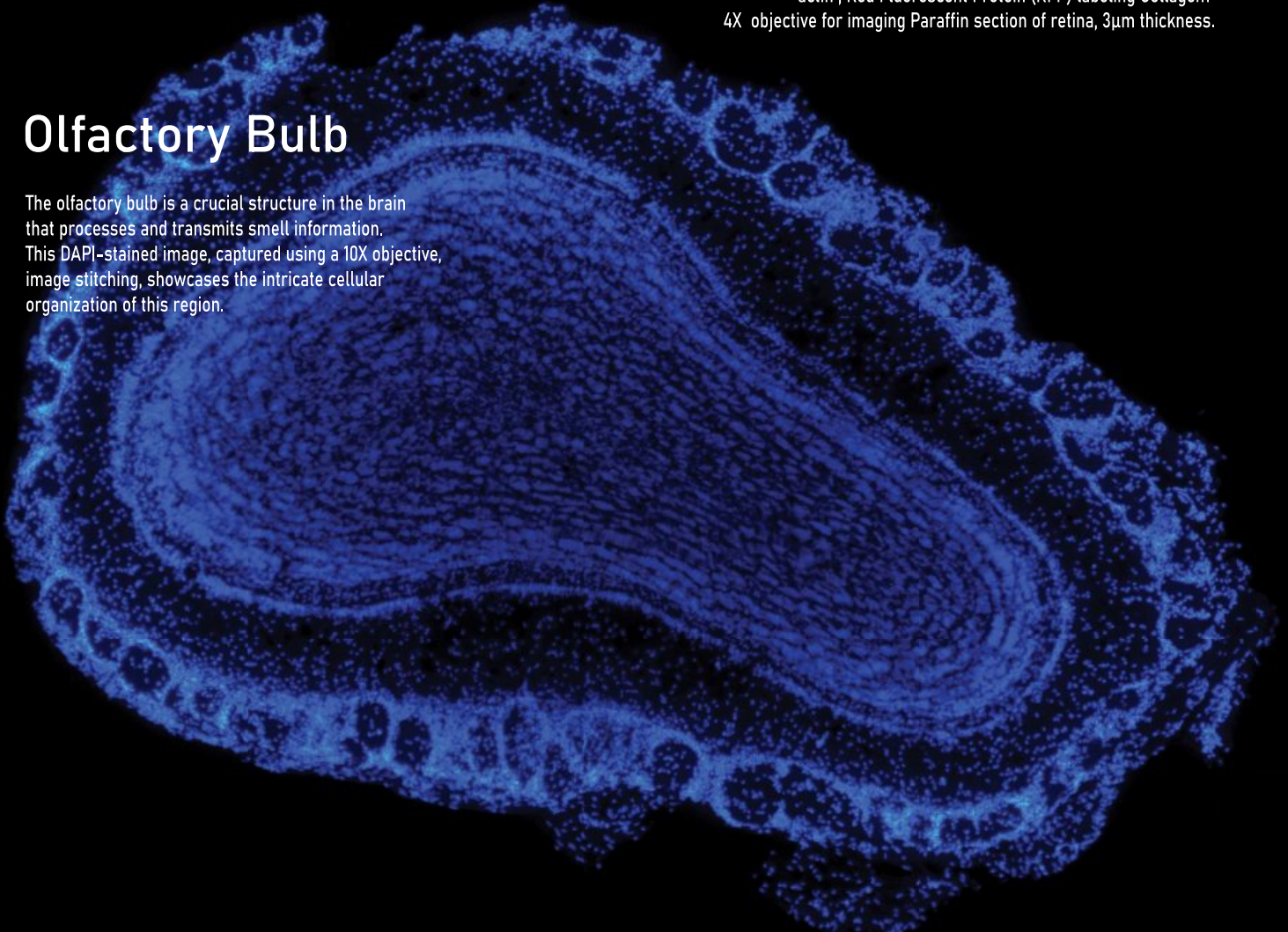

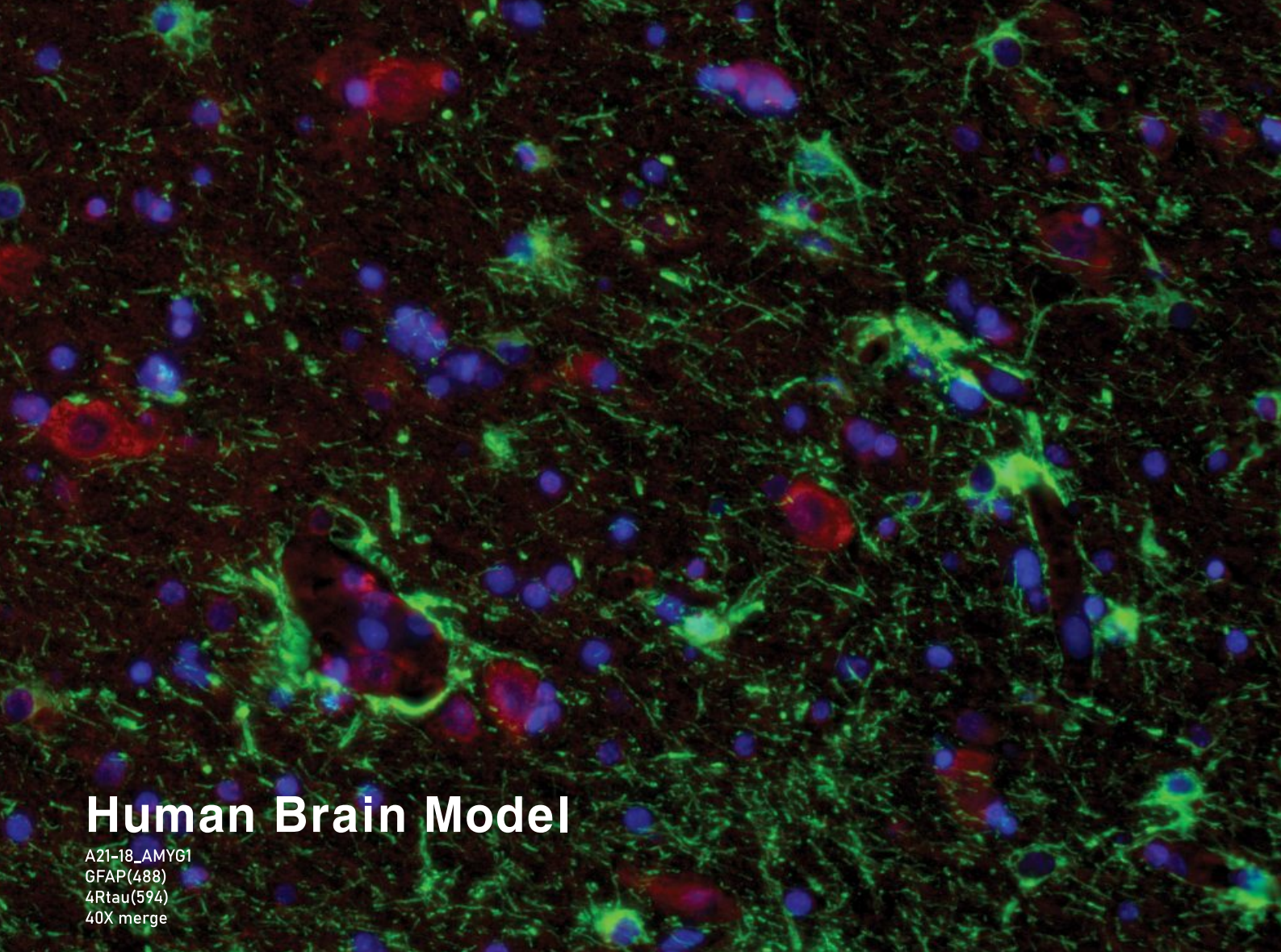


## Human Brain Model

A21-18\_AMYG1  
GFAP(488)  
4Rtau(594)  
40X merge

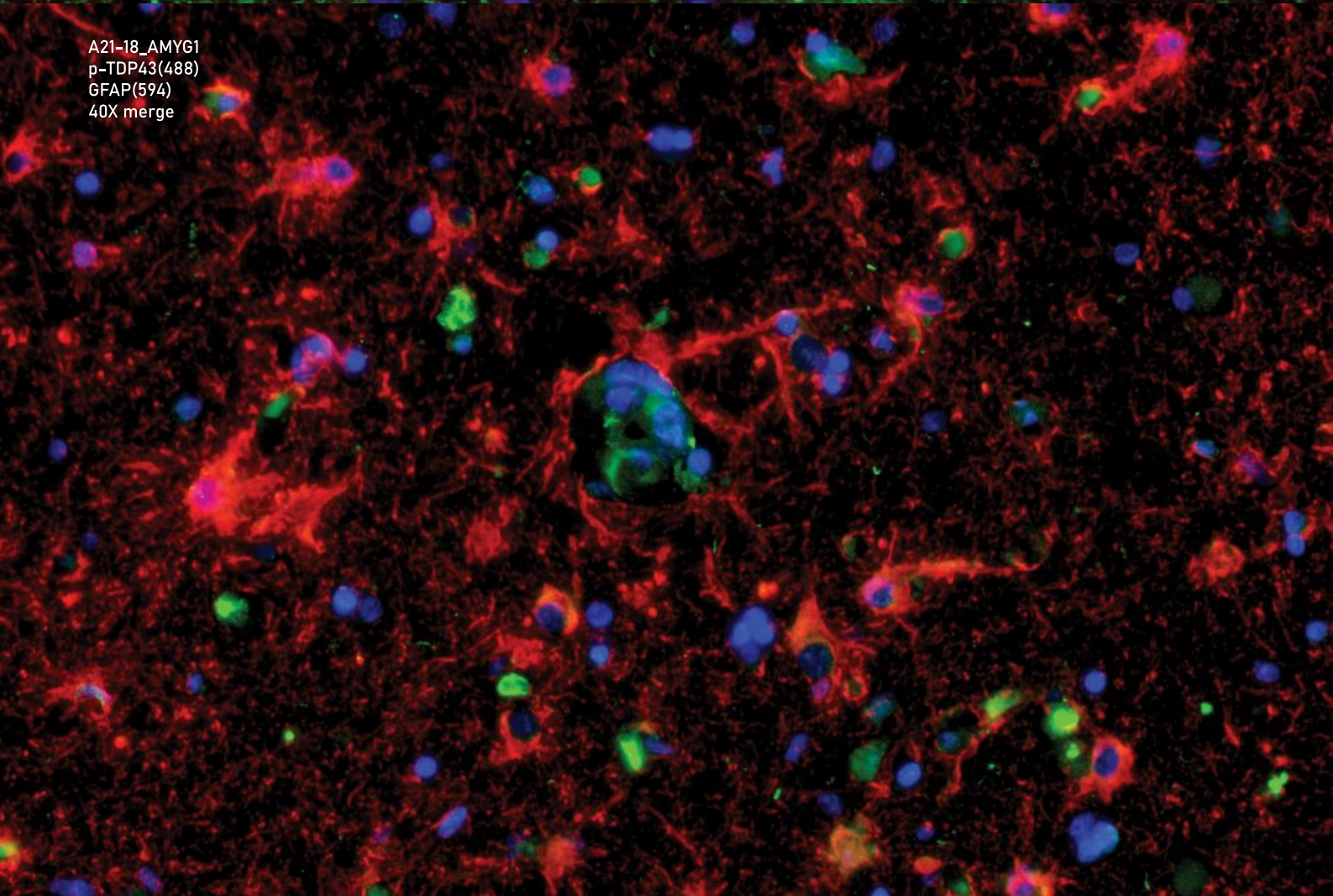


A21-18\_AMYG1  
p-TDP43(488)  
GFAP(594)  
40X merge

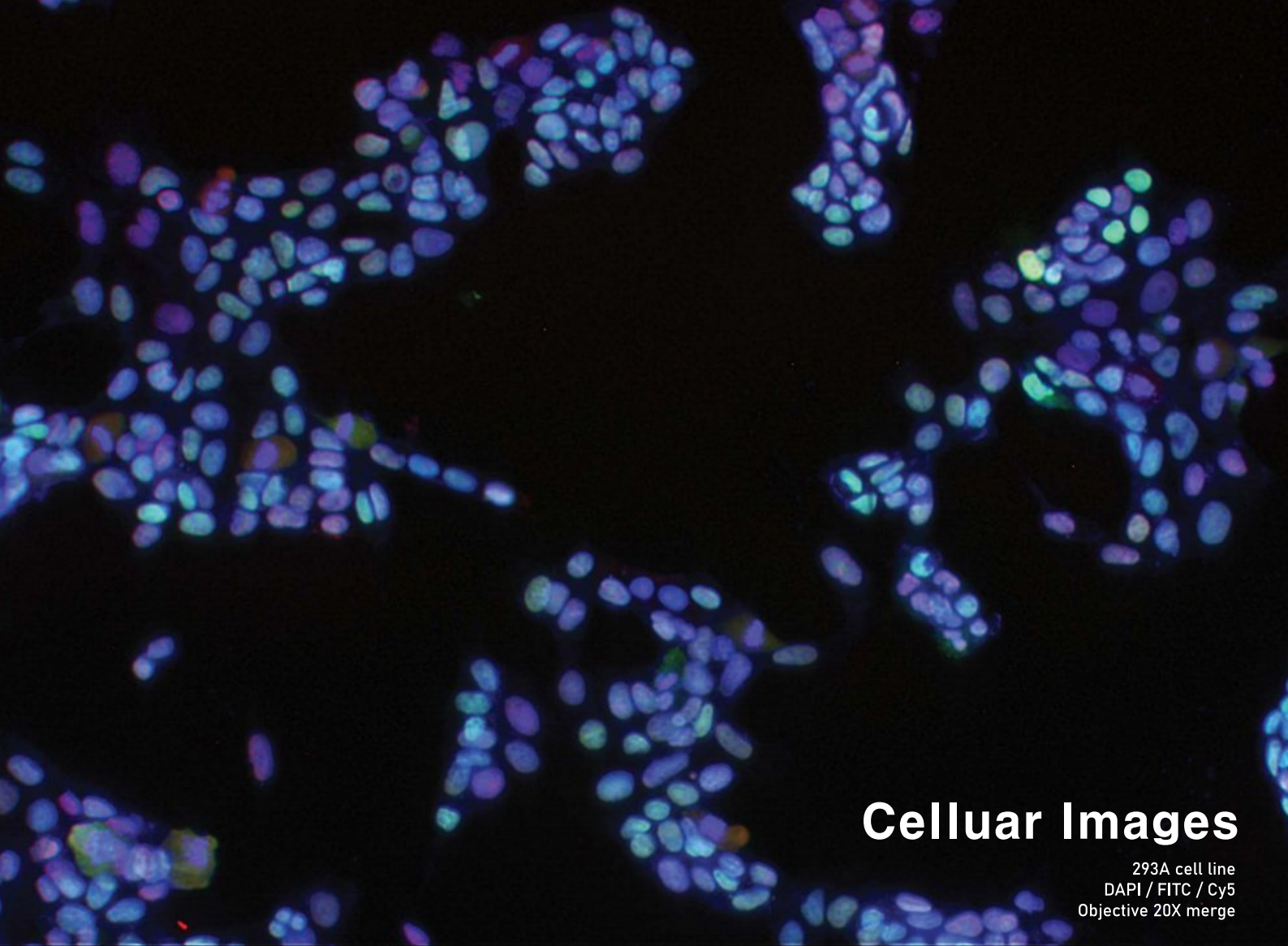


## Celluar Images

293A cell line  
DAPI / FITC / Cy5  
Objective 20X merge

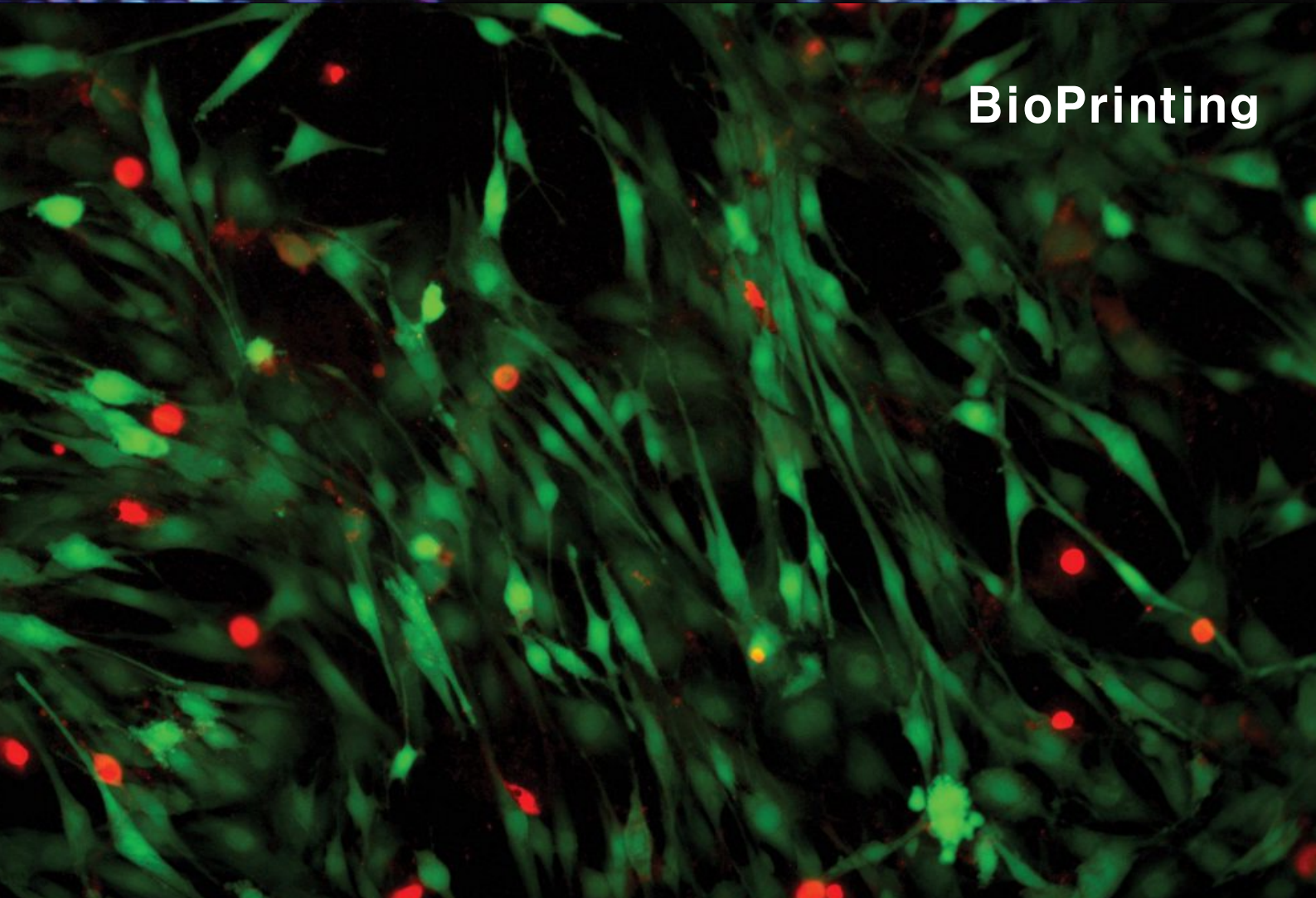



## BioPrinting

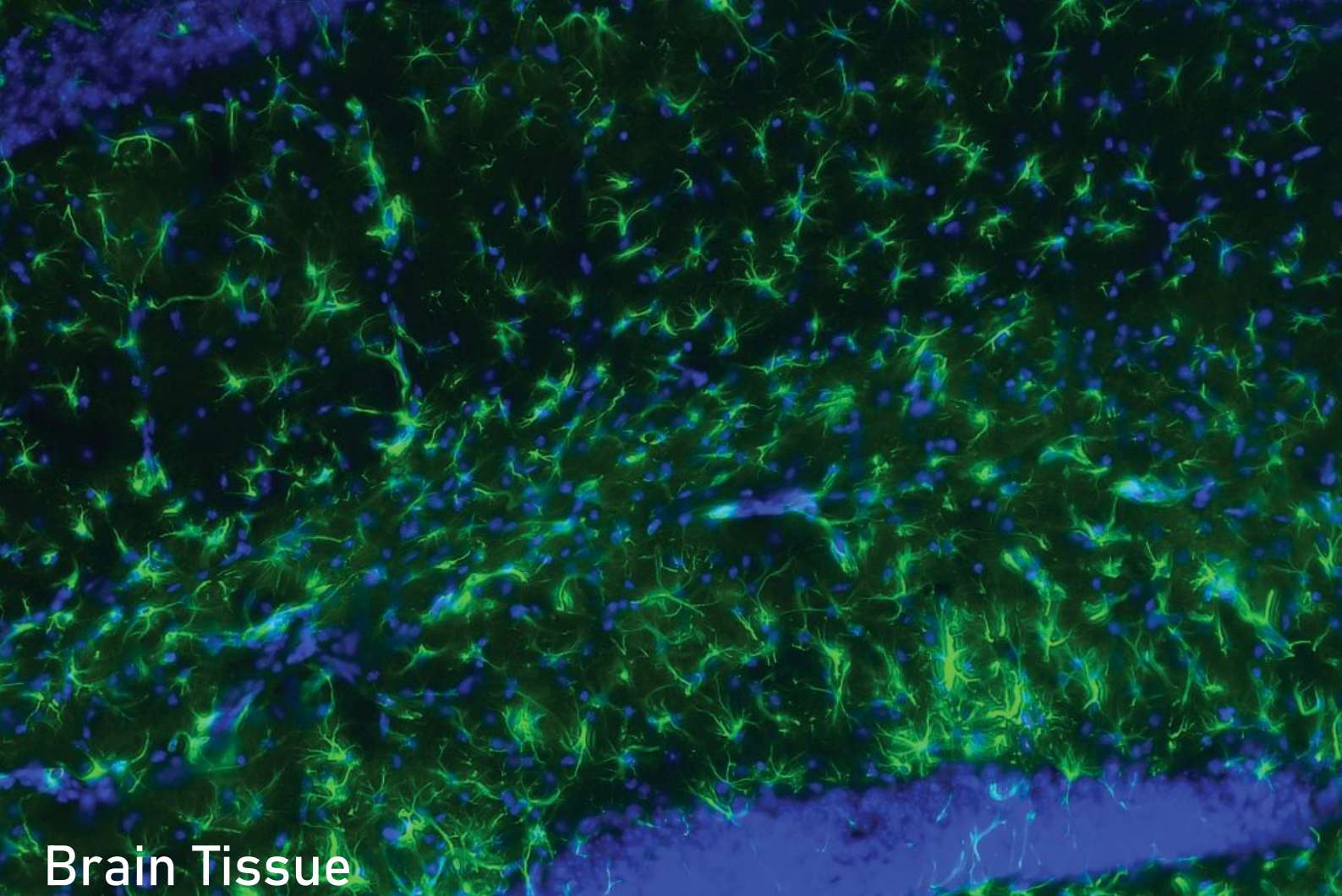

## Brain Tissue


Mouse brain tissue fluorescently stained for GFAP, photographed at 200x magnification. Fluorescent staining with DAPI, Alexa Fluor 488, and Alexa Fluor 568 was performed for neuroscience research

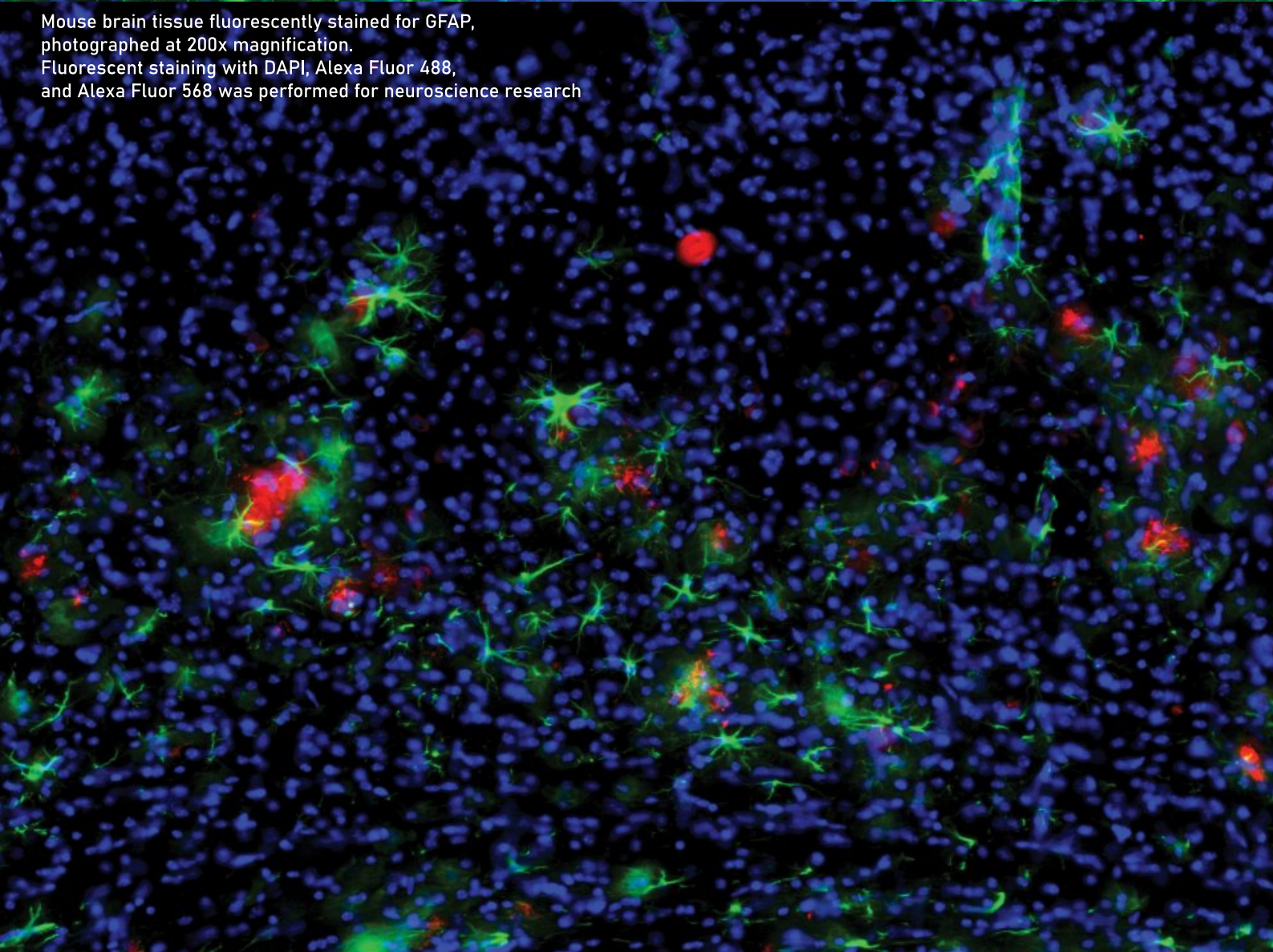

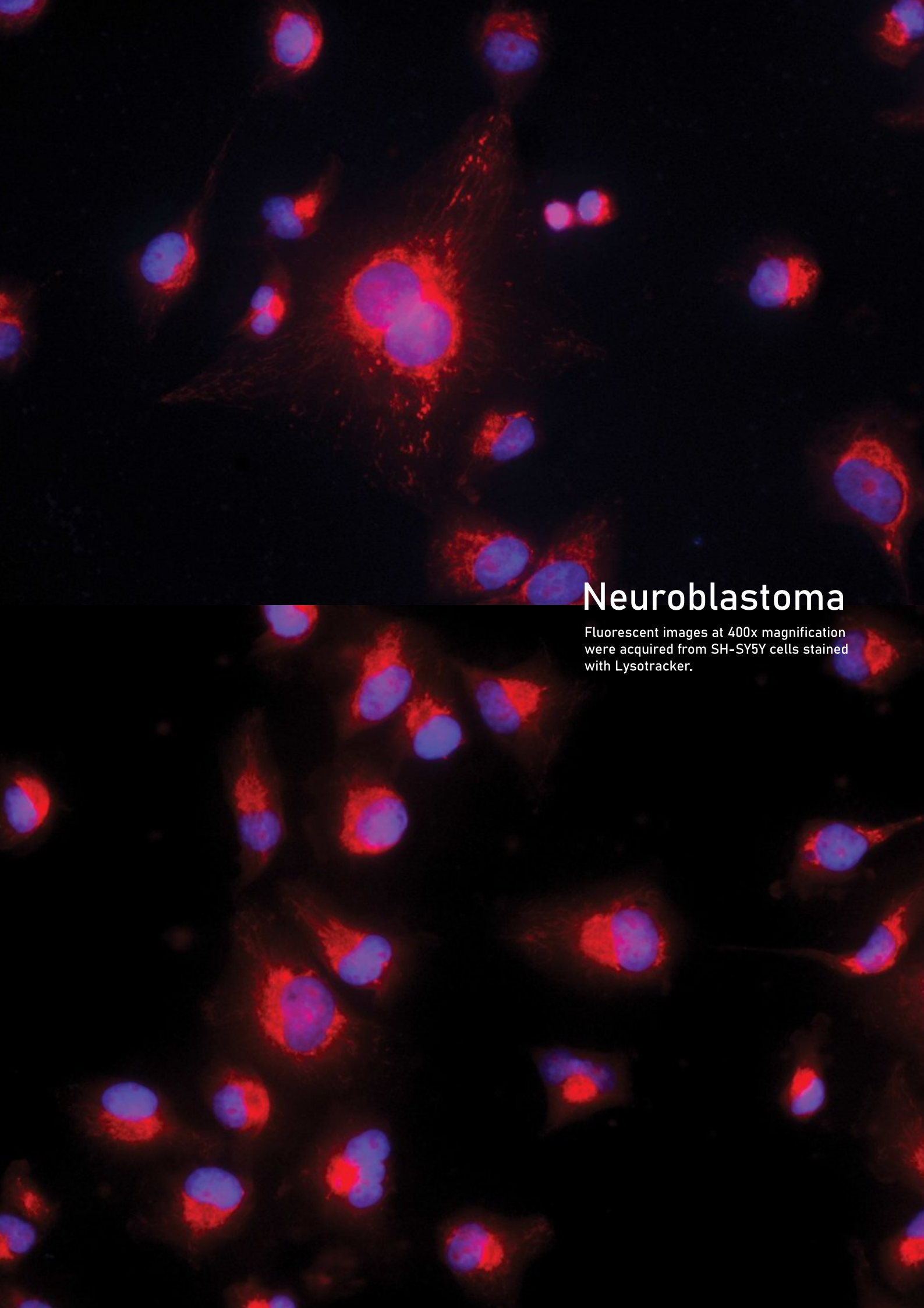

# Neuroblastoma

Fluorescent images at 400x magnification were acquired from SH-SY5Y cells stained with Lysotracker.

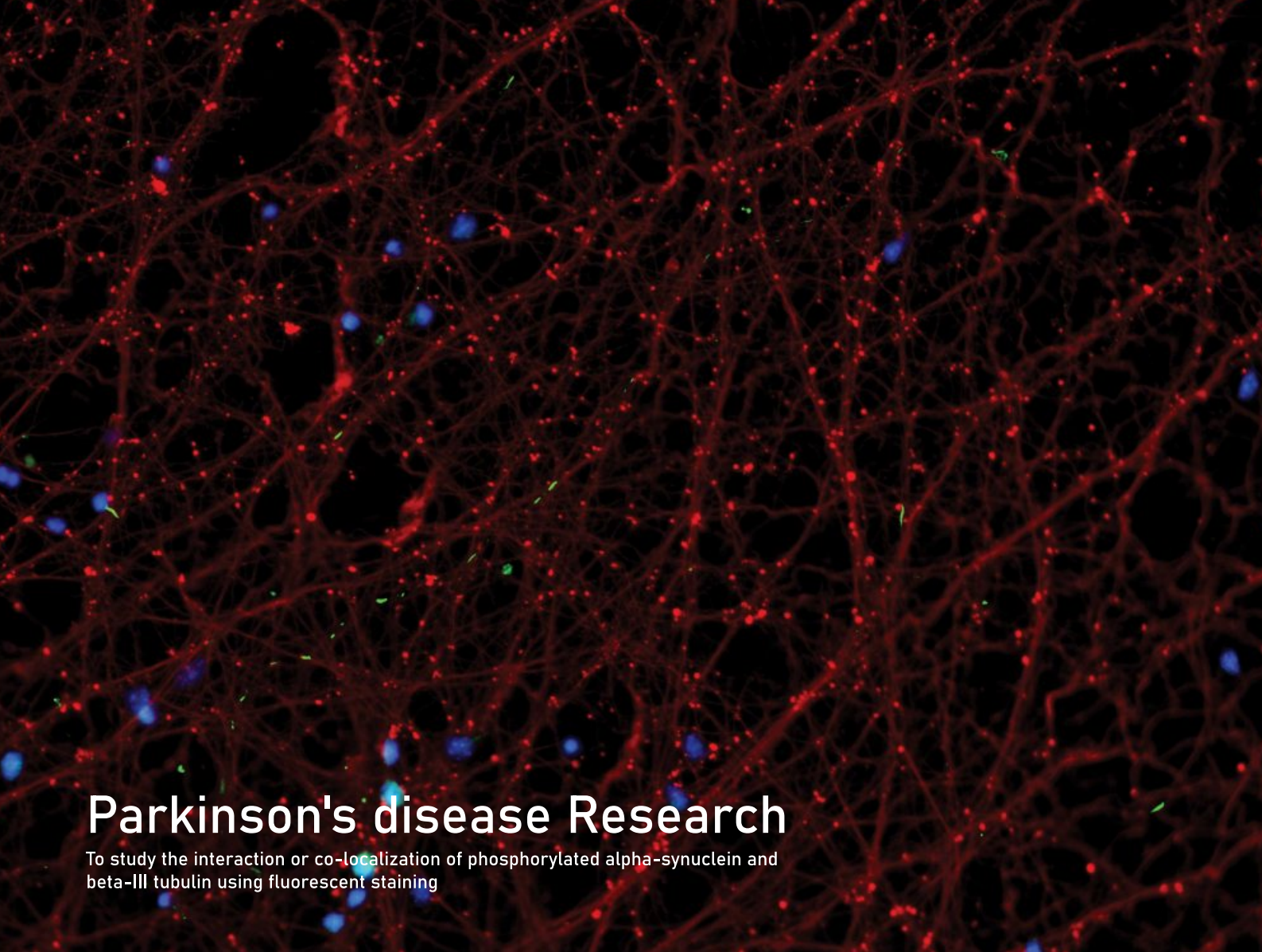



## Parkinson's disease Research

To study the interaction or co-localization of phosphorylated alpha-synuclein and beta-III tubulin using fluorescent staining

## Neurosphere


Neurospheres from primary cortical culture were fluorescently stained and imaged at 400x magnification


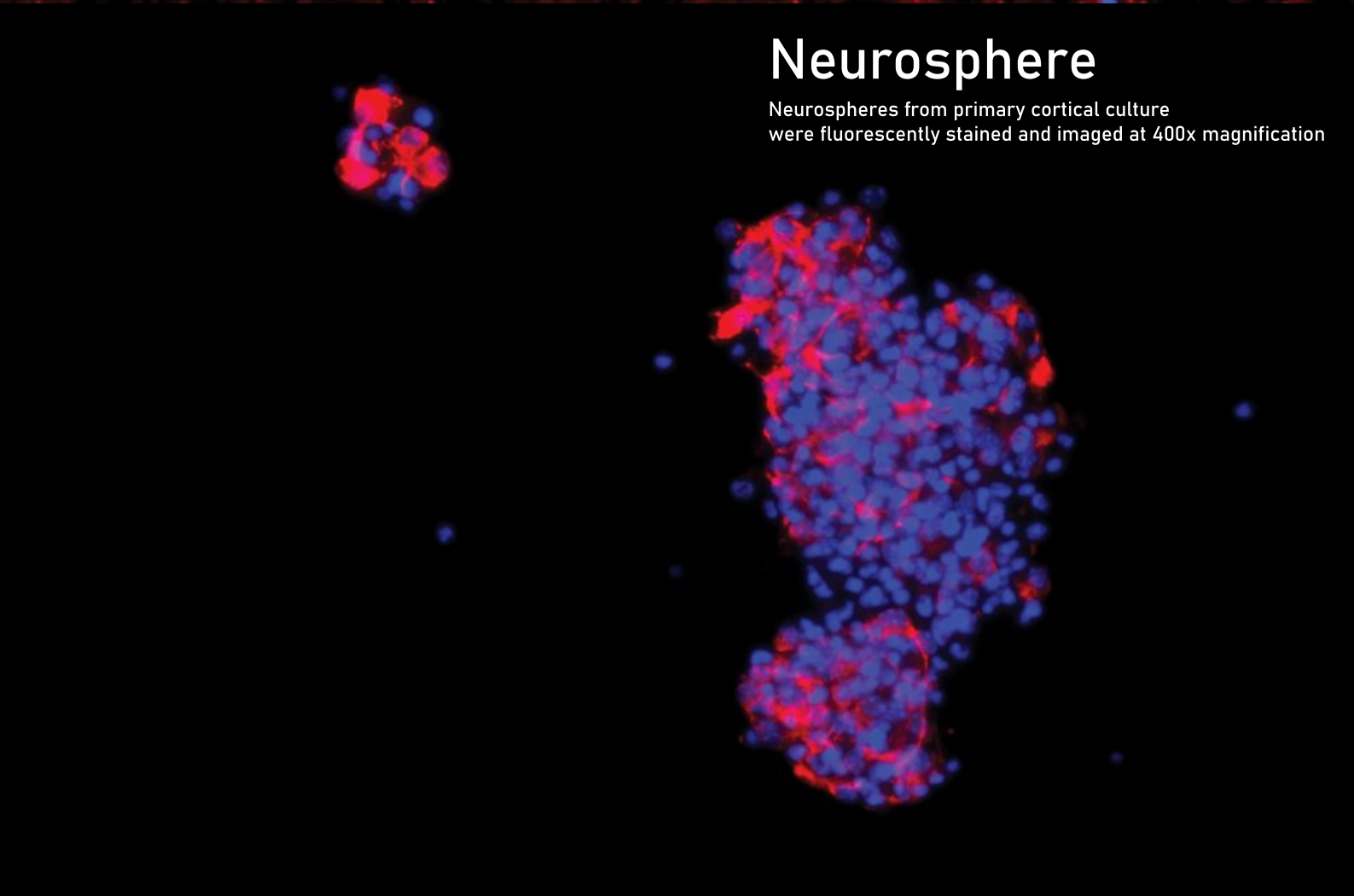


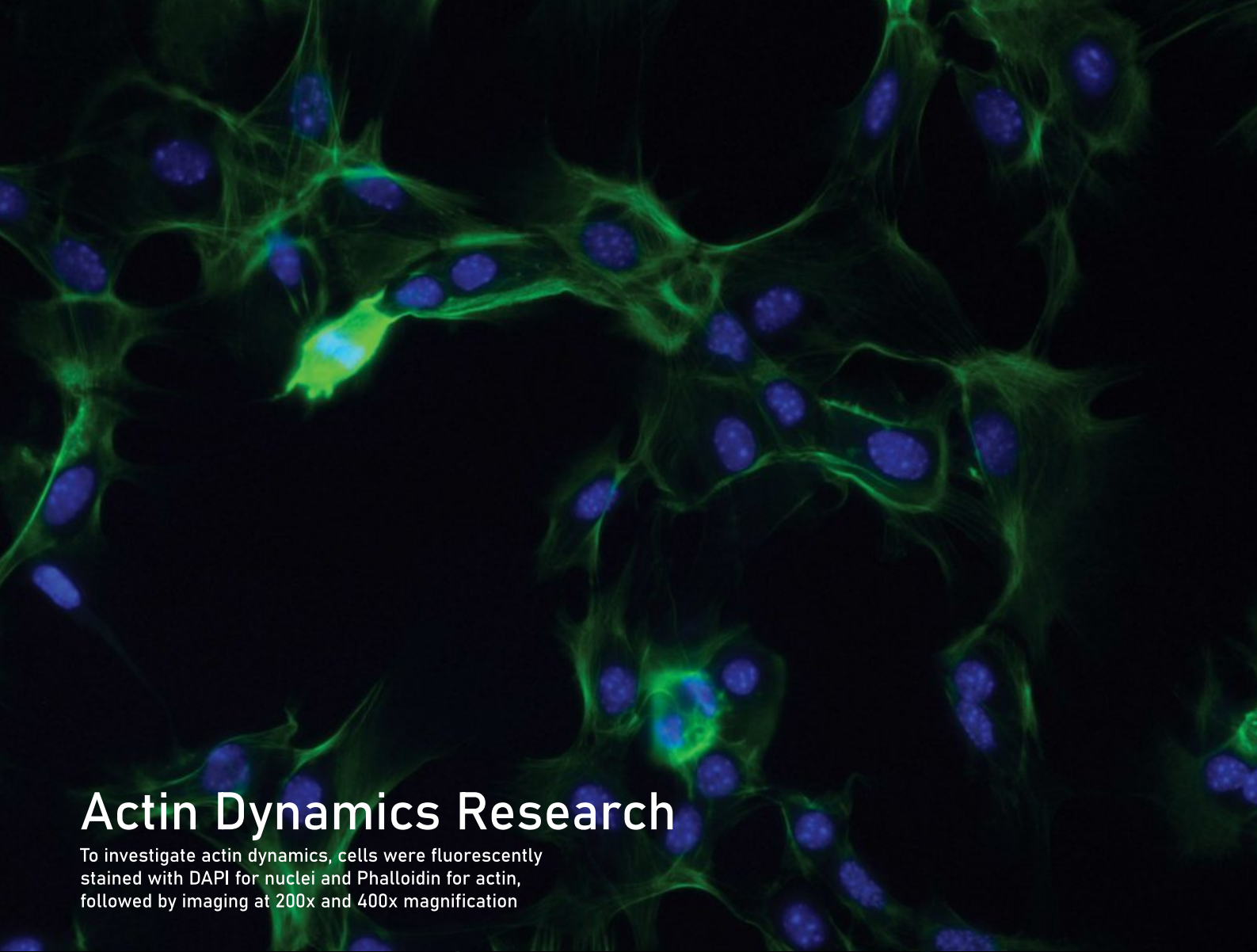

# Actin Dynamics Research

To investigate actin dynamics, cells were fluorescently stained with DAPI for nuclei and Phalloidin for actin, followed by imaging at 200x and 400x magnification

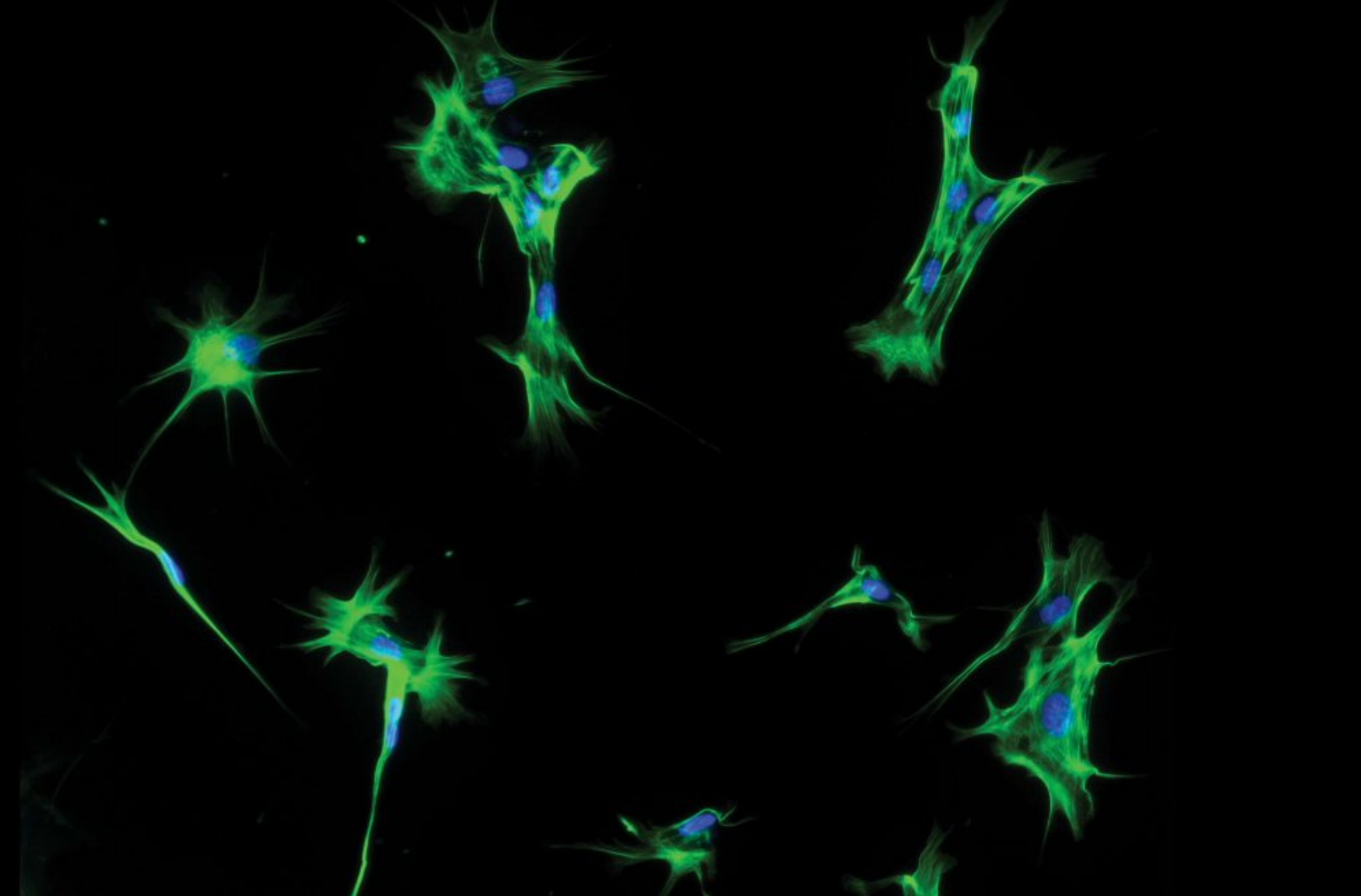

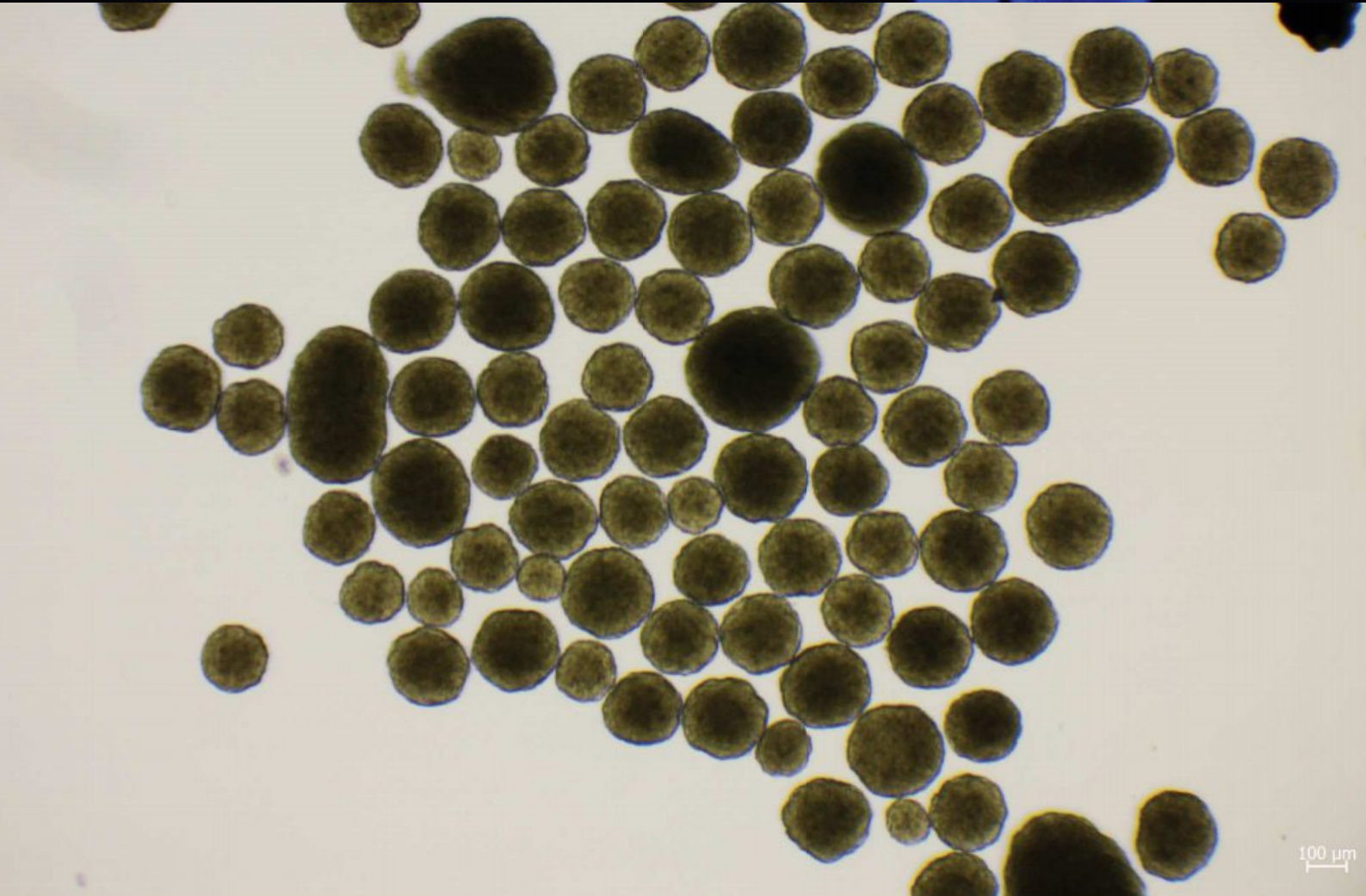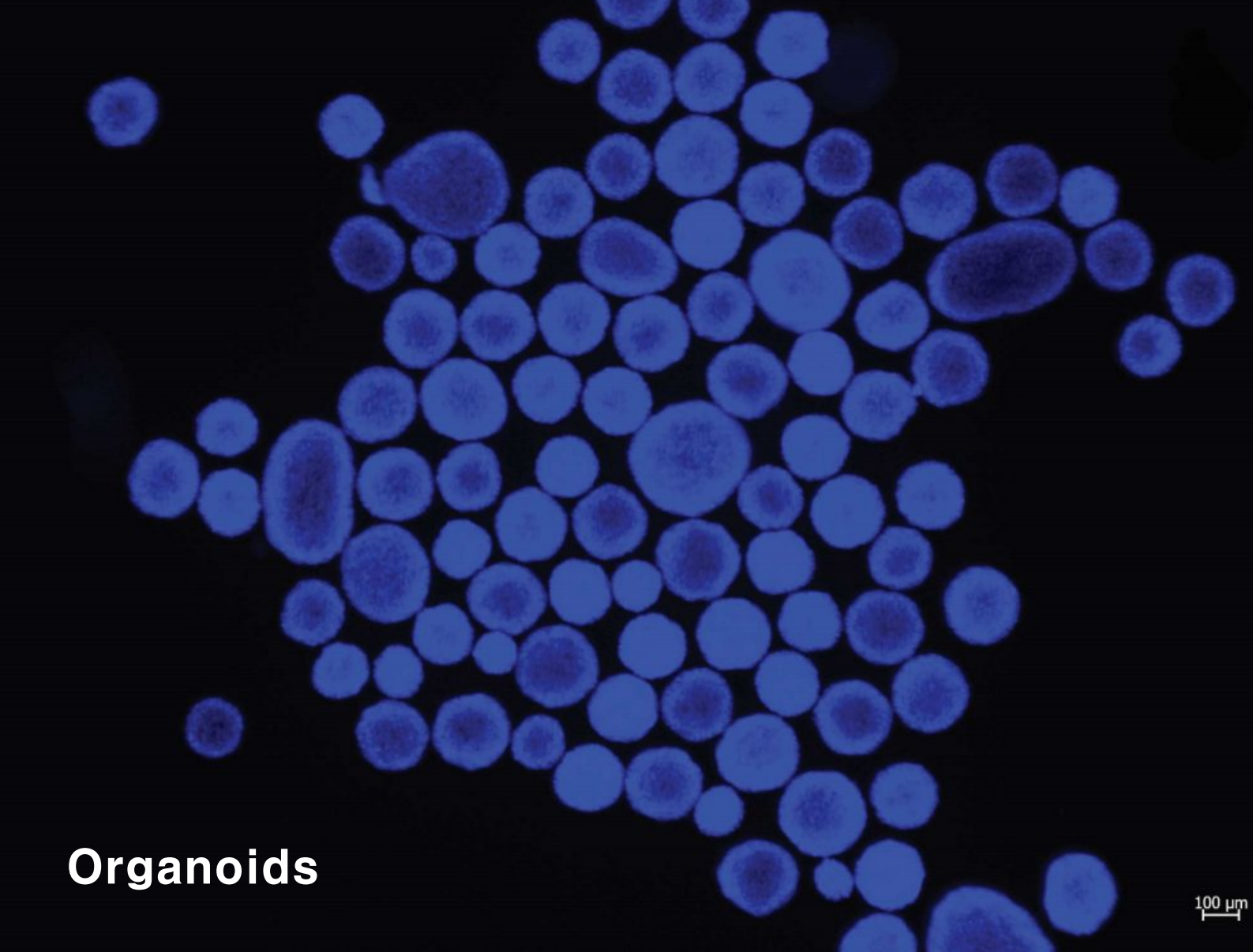




# Spheroids

TPBC(Triple-positive breast cancer) model / E-cadherin / Diameter 300um / 20X Z-Stack

Confocal microscope  
(JAPAN)

*Red*

Fluorescence microscope  
(Germany)

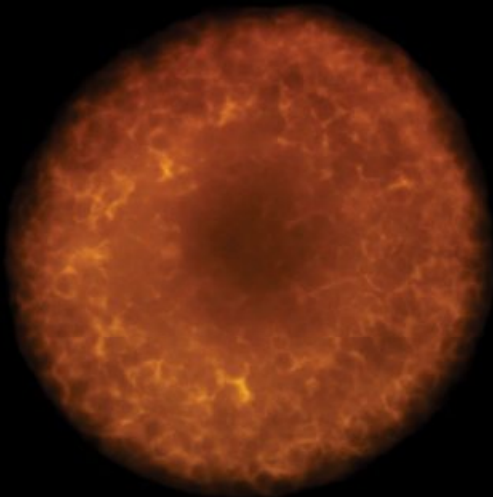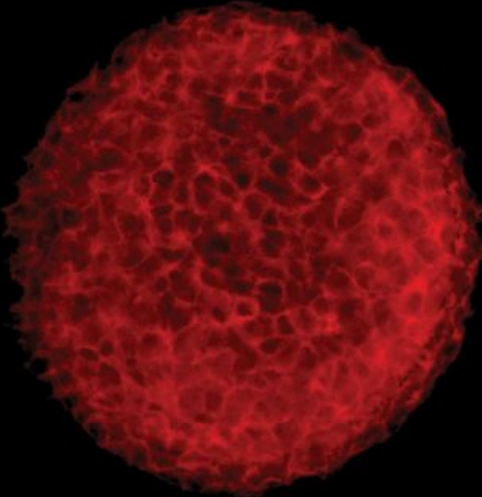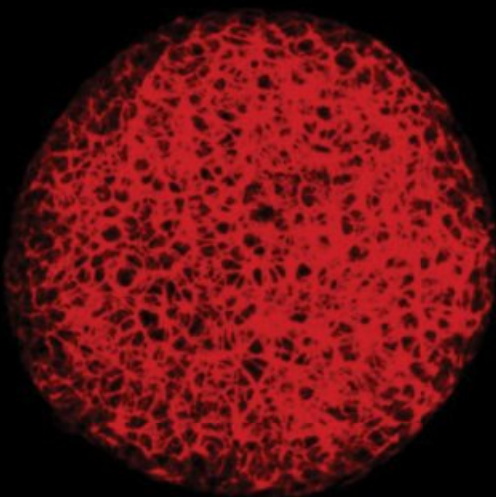





DAPI

Ki-67





E-cadherin

Merge




Confocal microscope  
(JAPAN)

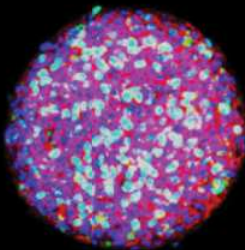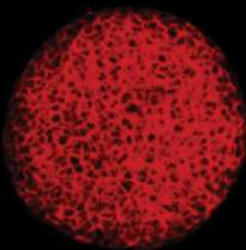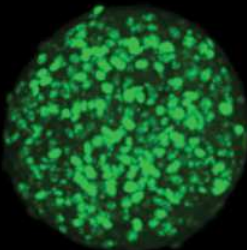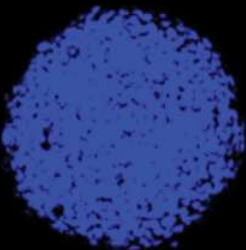

*Red*

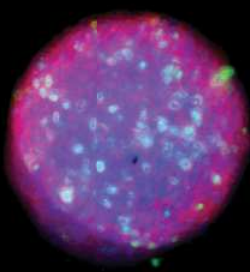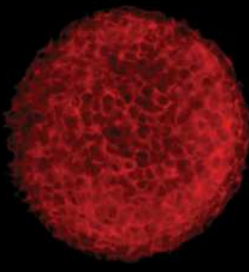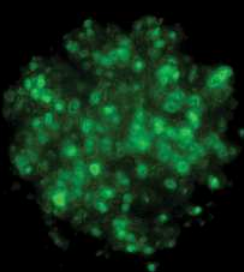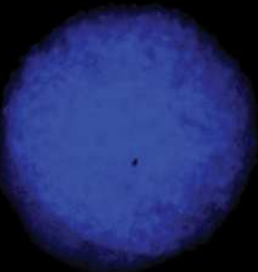

Fluorescence microscope  
(Germany)

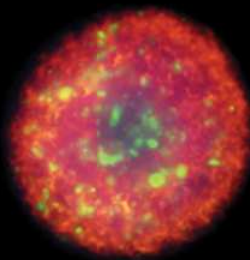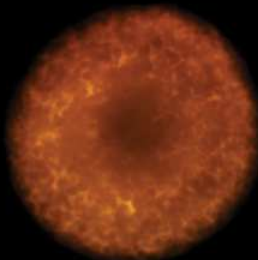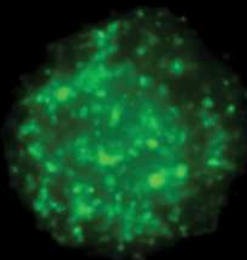


300  $\mu$ m

# Fluorescence Microscopy Fundamentals



## Easy-to-Operate Design

The Red microscope uses LED illumination that eliminates the need for bulb centering, providing a compact and user-friendly experience.

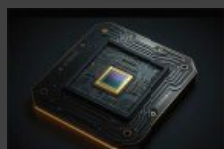

## High-Resolution Imaging

With a 12MP CMOS camera and objective lenses ranging from 1.25X to 100X, the Red delivers high-quality, detailed images.



## Comprehensive Filtering

A wide range of fluorescence filters, including Cy5, Texas Red, mCherry, Cy3, TRITC, GFP, FITC and DAPI are available for versatile imaging.

## Specification of *Red*


|                            |                         |                                                                                                                                                                                                                                                                                                                                                                               |            |            |         |
|----------------------------|-------------------------|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|------------|------------|---------|
| Microscope<br>Model : RF40 | Observation Method      | Bright field, Fluorescence, Phase contrast                                                                                                                                                                                                                                                                                                                                    |            |            |         |
|                            | Sample Holder           | Culture dish, Culture flask, 96well, 48well, 12well, 6well, plate, slide glase, confocal dish                                                                                                                                                                                                                                                                                 |            |            |         |
|                            | Objectives              | Plan fluor 4X, 10X, 20X, 40X (optional 1.25X, 2X, 60X) with 5-nospiece position turret                                                                                                                                                                                                                                                                                        |            |            |         |
|                            | Transmitted illuminator | Bulit-in transmitted white LED source (adjustable) with ND filter, Phase plate wheel                                                                                                                                                                                                                                                                                          |            |            |         |
|                            | Stage                   | Stage size: 277.5 x 210mm / X-Y mechanical stage moving range: 120x80mm                                                                                                                                                                                                                                                                                                       |            |            |         |
|                            | Focus                   | Focus knob resolution : 1 um                                                                                                                                                                                                                                                                                                                                                  |            |            |         |
|                            | Fluorescence parts      | High power 4 channel peak LED (wave length 365nm, 465nm, 550nm, 620nm )                                                                                                                                                                                                                                                                                                       |            |            |         |
|                            |                         | DAPI                                                                                                                                                                                                                                                                                                                                                                          | EX375/28nm | EM460/50nm | DM415nm |
|                            |                         | FITC                                                                                                                                                                                                                                                                                                                                                                          | EX480/30nm | EM535/40nm | DM505nm |
|                            |                         | Texas Red                                                                                                                                                                                                                                                                                                                                                                     | EX560/40nm | EM635/60nm | DM600nm |
|                            |                         | Cy5                                                                                                                                                                                                                                                                                                                                                                           | EX620/60nm | EX700/75nm | DM660nm |
| Camera                     | Sensor                  | 12 Megapixel, 1/1.8 inch, Sony IMX226 color sensor with USB3.0 interface                                                                                                                                                                                                                                                                                                      |            |            |         |
|                            | Resolution              | Maximum 4000x3000 / 2000x1500                                                                                                                                                                                                                                                                                                                                                 |            |            |         |
| Software                   | Software                | Red Image Analysis Software (Made by LE.AM Solution Inc.)                                                                                                                                                                                                                                                                                                                     |            |            |         |
|                            | Functions               | Capture, Save, Quick & Batch Save, BF mode, Fluorescence mode, R-G-B Manual & Auto Merge, Back ground black balance. Scale bar on the image, Fluorescence cell count, Auto Image intensity > CX/ CY/ Width / Height / Area / Intensity Min / Intensity Avg / Intensity Max. Video recording, Time lapse, Z-stack multiple images [ Image file Formats : TIFF, BMP, PNG, JPG ] |            |            |         |
| Environment                | Dimension               | W302 X D311 X H329 mm                                                                                                                                                                                                                                                                                                                                                         |            |            |         |
|                            | Weight                  | Approximately 25.6kg                                                                                                                                                                                                                                                                                                                                                          |            |            |         |
|                            | Power consumption       | 230.1 W                                                                                                                                                                                                                                                                                                                                                                       |            |            |         |
|                            | Electrocal Ratings      | Power Adaptor : INPUT AC 100-240V~, 50/60Hz, 3.5A / OUTPUT DC 19.5Vdc, 11.8A                                                                                                                                                                                                                                                                                                  |            |            |         |
|                            |                         | RF40 Equipment: INPUT DC in 19.5V, 11.8A                                                                                                                                                                                                                                                                                                                                      |            |            |         |

*Red*

Strong Fluorescence Imager

Sometimes,  
you don't need a  
confocal microscope



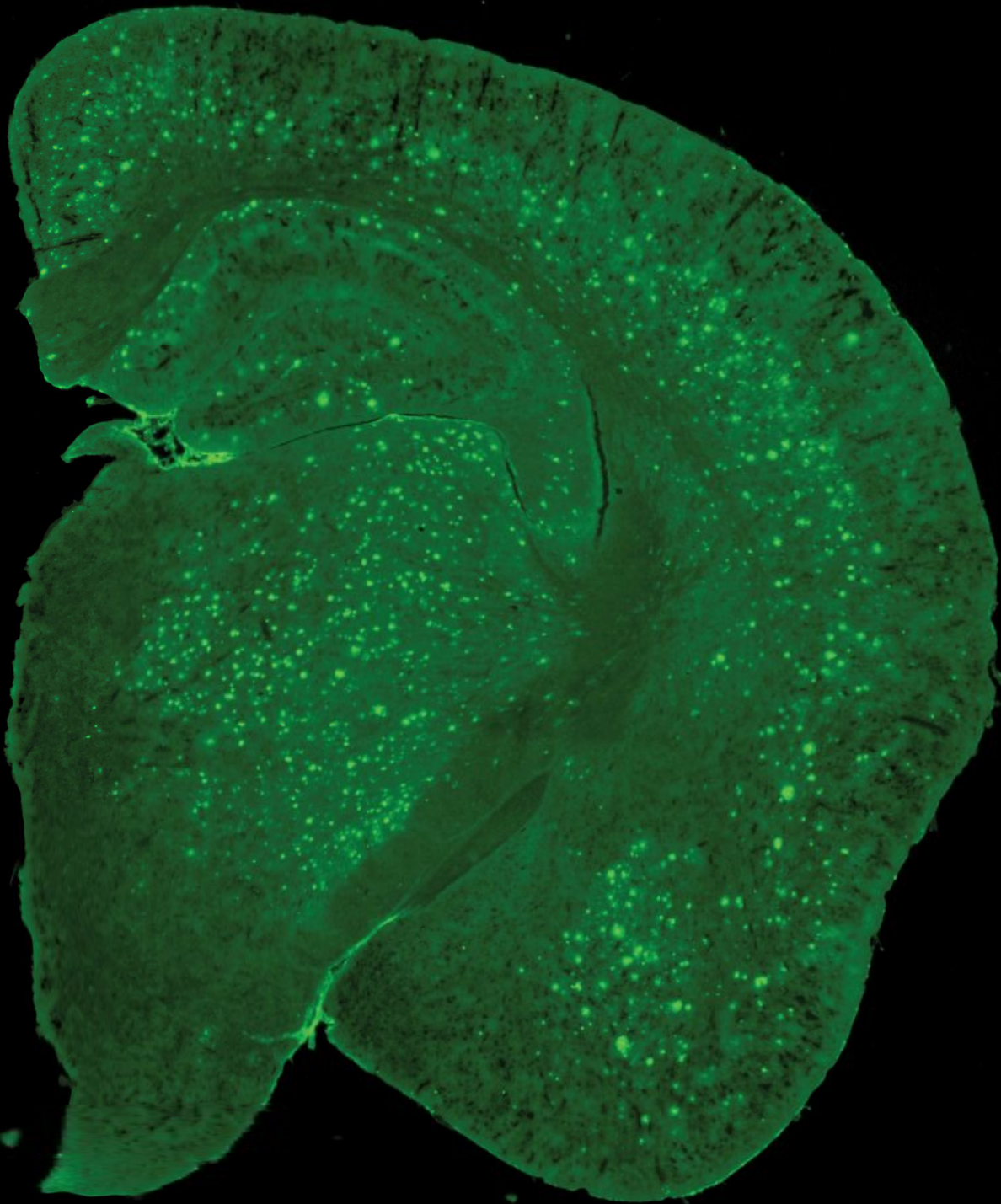

Made by **LE.AM** KOREA

LE.AM Solution Inc

Office : +82-31-434-2710

e-mail : leamsol@leamsol.com

www.leamsol.com

